UVL-4F Longwave UV Lamp

User's Manual

Important: Please read the instructions carefully prior to using the lamp in order to avoid malfunction or damage.

Operation

Easily switch between the longwave UV and the flashlight by using the selector located on the side of the lamp. Flashlight Off Longwave UV UV Tube Battery compartment

Insert 4 AA batteries (UM-3) not included with the lamp. To insert batteries, slide the battery cover out and flip upwards. Position the batteries as shown on the inside of the battery compartment. Slide back the battery cover.

Keep the lamp away from extreme temperature changes, excessive humidity, moisture, rain or other harsh environments. Do not mix old and new batteries together. It is not recommended to use the fluorescent lamp function when the energy level is low, it will blacken the fluorescent tube. Remove the batteries if the lamp is not to be used for a long period.




Replacement Parts Replacement Tube, 4W longwave BLB

Part Number 34-0010-01

Bulb cover

End cap

UV tube

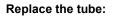


- 1. Remove the batteries prior to replacing the tube.
- 2. Pull off the end cap and circular bulb holder on the end of the unit, followed by the clear cover side panel.
- 3. Lightly twist the tube by the base to release the tube.
- 4. Replace with a 4 watt longwave tube.
- 5. Replace the clear cover.
- 6. Reconnect the end cap and bulb cover back into position.

Replacing the batteries:

When the tube becomes dim, change the batteries to retain lamp performance.

- 1. Slide the battery cover out and lift it up.
- 2. Remove the old batteries and replace with four new "AA" size "UM3" batteries, noting the battery orientation for proper insertion.
- 3. Slide the battery cover back into position.

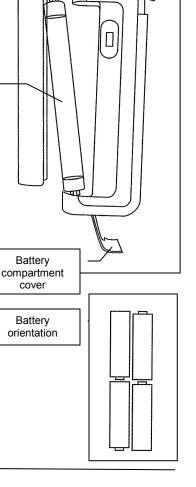


UVP, LLC 2066 W. 11th Street, Upland, CA 91786 Tel: (800) 452-6788 or (909) 946-3197 Fax: (909) 946-3597 Ultra-Violet Products Ltd. Unit 1, Trinity Hall Farm Estate, Nuffield Road, Cambridge CB4 1TG UK Tel: +44(0)1223-420022 Fax: +44(0)1223-420561 Web Site: www.uvp.com

UVL-4F Longwave UV Lamp

IsunsM s'192U

using the lamp in order to avoid malfunction or damage. Important: Please read the instructions carefully prior to


Operation

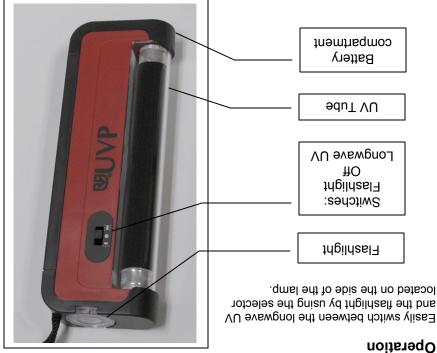

inside of the battery compartment. Slide back the battery cover. the battery cover out and flip upwards. Position the batteries as shown on the Insert 4 AA batteries (UM-3) not included with the lamp. To insert batteries, slide

the lamp is not to be used for a long period. energy level is low, it will blacken the fluorescent tube. Remove the batteries if together. It is not recommended to use the fluorescent lamp function when the moisture, rain or other harsh environments. Do not mix old and new batteries Keep the lamp away from extreme temperature changes, excessive humidity,

Euq csb Remove the batteries prior to ۱. Replace the tube: Bulb cover Replacement Tube, 4W longwave BLB Part Number Replacement Parts

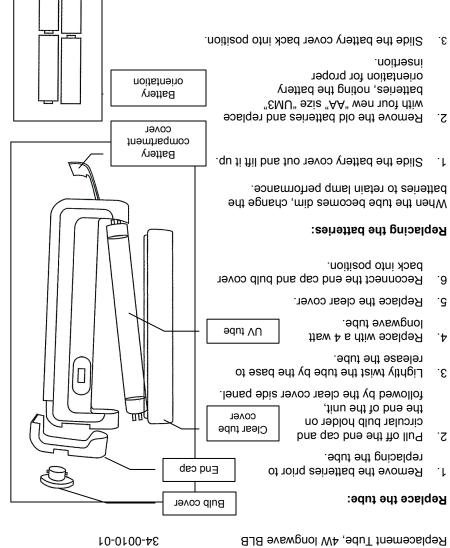

moo.qvu.www :9ji2 d9W Cambridge CB4 1TG UK Tel: +44(0)1223-420022 Fax: +44(0)1223-420561 Ultra-Violet Products Ltd. Unit 1, Trinity Hall Farm Estate, Nuffield Road, 7925-349 (909) :xb7 7915-349 (909) 946-3597 **UVP, LLC** 2066 W. 11th Street, Upland, CA 91786 Tel: (800) 452-6788 or

3.

.2

.9

.6

.4